BRESSER®

- DE** Bedienungsanleitung
- GB** Instruction Manual
- FR** Mode d'emploi

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General Information

About this Instruction Manual

Please read the safety instructions in this manual carefully. To avoid damage to the unit and the risk of injury, only use this product as described in the manual. Keep the instruction manual handy so that you can easily look up information on all the functions.

DANGER!

You will find this symbol in front of every section of text which deals with the risk of severe injury or even death in the event of improper use.

CAREFUL!

You will find this symbol before every section of text that deals with the risk of minor to severe injuries resulting from improper use.

NOTE!

You will find this symbol in front of every section of text which deals with the risk of damage to property or the environment.

Intended Use

This product is intended only for private use. It was developed for the magnified display of things in nature.

General Warning

DANGER!

Tools with sharp edges and points are often used when working with this device. For this reason, store this device and all accessories and tools in a location that is out of the reach of children. There is a RISK OF INJURY!

DANGER!

Do not expose the device to high temperatures. Use only the supplied power adapter. Do not short circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short circuit, fires or even explosions.

DANGER!

Never bend, pinch, pull or run over the power and connecting cables or extensions/adapters. Protect the cables from sharp edges and heat. Before operation, check the device, cables and connections for damage.

Never use a damaged unit or a unit with damaged power cables, etc. Damaged parts must be exchanged by an authorized service centre immediately.

**DANGER!**

Children should only use the device under adult supervision. Keep packaging material (plastic bags, rubber bands, etc.) out of the reach of children! There is a RISK OF CHOKING!

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after use. In the event of contact with the eyes or mouth, rinse thoroughly with water. In the event of pain, contact a doctor immediately and take the substances with you.

**DANGER!**

Children should only use the device under adult supervision. Keep packaging material (plastic bags, rubber bands, etc.) out of the reach of children! There is a RISK OF CHOKING!

**CAREFUL!**

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after use. In the event of contact with the eyes or mouth, rinse thoroughly with water. In the event of pain, contact a doctor immediately and take the substances with you.

**NOTE!**

Do not disassemble the device. In the event of a defect, please contact your dealer. He will contact the Service Centre and can send the device in to be repaired, if necessary.

Do not expose the device to temperatures above 40°C.

Position your device so that it can be disconnected from the power supply at any time. The wall socket should always be located near the device and be easily accessible, since the plug on the power cord serves as a disconnecting device for the power supply.

All parts:

- ① Eyepieces
- ② Diopter adjustment ring
- ③ Binocular viewing
- ④ Objective
- ⑤ Objective clamping screw
- ⑥ Focussing wheel
- ⑦ Illumination
- ⑧ Guidance of microscope head
- ⑨ Power switch
- ⑩ Microscope stage with object plate (black/white)
- ⑪ Object clips
- ⑫ Fuse holder (with fuse)
- ⑬ Power cord
- ⑭ Fastening screw for guard cover
- ⑮ Guard cover

1. General/Location

Before starting to set up your microscope, select a suitable location for it. First make sure that your microscope is put up on a stable, vibration-free basis. In order to observe using the electrical illumination, a mains socket (230 V, 50 Hz AC) is needed.

2. Illumination

For the start-up, connect the power cord (13) with the mains socket. Turn on the illumination with the power switch (9) at the top of the microscope.

3. Observation

Place the object to be viewed in the centre of the microscope stage (10). If it is possible, clamp it on the edges with the object clips (11).

Now, look through the eyepieces of the binocular viewing (3) and change the spacing between the eyepieces (1) so that you get a round, clear field of view. The definition is adjusted by slowly turning the focussing wheel (6).

If you have to observe a very large or very small object, you can adjust the microscope. Loosen the holding screw on the back and move appropriately the microscope head on its guidance (8). If you have found the correct position, tighten the holding screw again.

4. Diopter adjustment

Turn the diopter adjustment ring (2) of the right eyepiece to the middle position (silver marking ring). Shut your right eye and look with the left eye through the left eyepiece. With the focussing wheel (6), focus correctly for your left eye. Now, look with your right eye through the right eyepiece. If the image isn't sharp, adjust the focus for your right eye by turning the diopter adjustment ring (2). Now, the microscope is correctly adjusted.

5. Exchange of the lamp

Make sure that the microscope is separated from the mains. Let the lamp cool down, if necessary, since lamps can become very hot in operation.

Loosen the screw (14) and drag the guard cover (15) out. Now, you can easily remove the lamp. Insert a new lamp (12 V, 10 W) in the lamp holder. Grasp the new lamp not with your fingers, do this best with a cloth. Now, you can back up the guard cover (15) and tighten the screw (14) again. Tighten the screw only in a hand-screwed way.

6. Exchange of the fuse

The microscope is equipped with two fuses (12). First make sure that the microscope is separated from the mains. To replace a fuse turn out the fuse holder using a small screwdriver. Remove the defective fuse and insert a new one (0.5 A, 250 V). Screw the fuse holder back into the device.

7. Care and maintenance

Cleaning the lenses: Sweep the lens with a lens tissue or soft fabric immersed with a small amount of alcohol (e.g. isopropanol).

Cleaning the painted parts: The dust on the painted parts can be removed by gauze. For the grease spots, the gauze immersed slightly with aviation gasoline is recommended. Do not use

organic solvents such as alcohol, ether or other thinner etc. for cleaning the painted parts or plastic components.

Your microscope is a high-quality optical instrument. Avoid disassembling the microscope. Improper disassembling affects the high utility of the microscope. Moreover, the warranty expires.

Cover the microscope with the dust protective cover and place it there, where it is dry and dust-free. We suggest the storage of all objectives and eyepieces in a closed container with drying agent.

8. Exchange of objectives (optional)

To change the objectives, loosen the objective clamping screw (5). Now, you can sideways pull out the objective. Then insert another objective. Mount the objective clamping screw (5) again and tighten it in a hand-screwed way, so slipping out of the objective is excluded.

9. Technical data

Binocular viewing

Eyepiece: WF 10x (Ø 30.5 mm)

Objective: 2x

Magnification: 20x

Illumination: 12 V, 10 W (bulb)

Fuse: 0.5 A, 250 V

Accessories: Spare lamp, spare fuse, wrench, 2 rubber eyecups, dust protecting cover, instruction manual.

10. Only for EU countries

Do not dispose of electric equipment together with household waste material!

In observance of European Directive 2002/96/EC on waste electrical and electronic equipment (WEEE) and its implementation in accordance with national law, electric equipment that have reached the end of its life must be collected separately and returned to an environmentally compatible recycling facility.

Discharged batteries and damaged re-chargeable batteries must be disposed of at special battery collection points. Information is available from your local disposal agent or local authority regarding the disposal of devices or batteries.

11. Declaration of Conformity

Bresser GmbH has formed a „Conformity Declaration“ in line with the applicable guidelines. This declaration is available for review.

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Un « Certificat de conformité » conforme aux directives applicables, établi par la société Bresser GmbH est disponible.